

J6-MI Centrifuge Users

- Contact Julie (jungmi@uoguelph.ca)
- Sign the Log book
- Balance the rotor
- Read the instruction
- Clean after use

NORMAL RUN MODE



For low-temperature work, refrigerate the rotor and sample before the run. POWER must be ON before the chamber door can be opened or closed. Carefully lower the rotor straight down over the drive spindle, tighten the knob, and close the chamber door.

TO START A MANUAL RUN:

1. Turn the key switch to NORMAL.
2. Press [ROTOR]. Press the rotor entry code for the rotor in use (see table below). Press [SPEED]. Enter the required speed. Press [TIME]. Continue and enter TIME, TEMP, and ACCEL, DECEL, and COMP ADJ if necessary.
3. Press [ENTER/RECALL].
4. Press [START].

RPM	JS-4.2 an JS-4.2A * g-force
500	71
1 000	284
1 500	640
2 000	1 140
2 500	1 780
3 000	2 560
3 200	2 910
3 500	3 480
4 000	4 550
4 200	5 020